

Myco 5-in-1 PLUS: Detect SIX mycotoxins with ONE extraction.

VICAM™

A Waters Business

1

Weigh 5 g of finely-ground sample into a extraction tube, then add 25ml of Aqua premix.



2

Extract the sample by vortexing at high speed for 2 minutes.



3

Filter the extract into a clean tube.



Afla-V™
QUANTITATIVE STRIP TESTS

Ochra-V™
QUANTITATIVE STRIP TESTS


Fumo-V™
QUANTITATIVE STRIP TESTS

Zearala-V™
QUANTITATIVE STRIP TESTS

T2/HT2-V™
QUANTITATIVE STRIP TESTS

DON-V™
QUANTITATIVE STRIP TESTS

Pipette 100 µL of extract to the sample well on the strip at 1 drop/second.

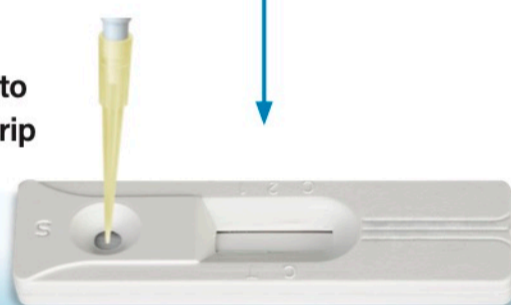

Place DON-V™ strip into the mini-incubator at 40 °C and close the lid. Incubate strip for 2-10 minutes. This can be started while doing the the sample extraction and filtration.



Pipette 100 µL of extract to the sample well on the DON-V strip at 1 drop/second and wait for the strip to develop for 5 minutes, leaving it inside the incubator (40 °C).



Solvent-free extraction!

Eliminate the expense and environmental hazards of toxic organic solvents.

Place the strip on a flat surface for 5 minutes.

While waiting for 5 min, prepare Vertu calibration by scanning the corresponding barcode for the specific mycotoxin kit.



Place the strip inside the Vertu reader (the round opening facing inwards), then press the center key.



To print result, move the cursor to 'P' using the right arrow key, then press the center key.

For more information visit www.vicam.com

VICAM, A Waters Business Headquarters:

34 Maple Street Milford, MA 01757, USA ■ Tel.: +1 800 338 4381 ■ +1 508 482 4935 ■ Fax.: +1 508 482 4972

Waters

THE SCIENCE OF WHAT'S POSSIBLE.™